


1. Use the standard Leica SP8 manual to operate the confocal
2. Before making use of the confocal, check the following elements on microscope stand:

- A. Lens must be placed inside the stand.
- B. Move switch to bright field.
- C. Height of condenser can be adjusted to increase quality of transmission image.

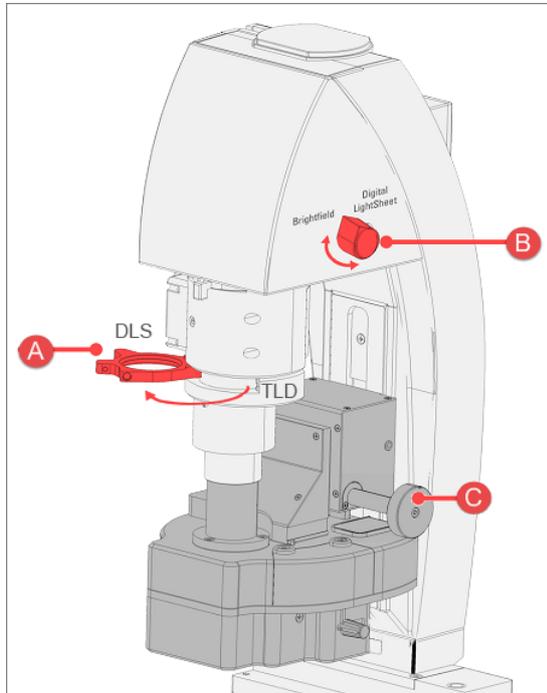

3. Select TLD option in Fluo Turret when obtaining transmission image.



CONTACT INFORMATION

For direct support at the microscope call **31105**.

For other question contact us via:

Martijn de Gruiter tel. 31105 h.degruiter@erasmusmc.nl