

Zeiss LSM900 Room Be311



1





A few points of attention:

- Do not use too much oil
- Do not use oil on the dry-lenses
- Clean the oil-lenses (40X, 63X) after use with a tissue and ethanol
- Do not eat or drink behind the microscope
- Throw away your garbage
- Be careful with the microscope

2



Turning on the system

- 1. Turn on the 'system'-switch behind the table, box above the box with the key. DO NOT TOUCH OTHER SWITCHES OR THE KEY!
- 2. Turn on the connector box on the microscope-table. DO NOT TURN ON OR OFF THE MICROSCOPE.
- 3. Turn on the computer
- 4. Double click 'Zen Blue'. If the whole system has been turned on click 'Yes' when asked to calibrate. The calibrate question will not appear if the system is on and only the software is started.

Looking and searching through the microscope

- 1. First use the 5X-, 10X- or 20X-lens!!
 - a. Choose a lens by clicking on the right lens in the 'Microscope'-window on the right of the screen.
- 2. In the Zen-program click 'locate'
- 3. On the microscope touch screen touch 'load position'
- 5. In the Zen-software click on one of the excitation wavelenghts (DAPI, 488 561). This will automatically choose the correct settings and open the shutter.
- 6. At 'Reflected light' you can close the shutter with 'Off'
- 7. Push the slider on the microscope in (right side of the oculairs) and focus your slide
- 8. If necessary change to the 40X- or 63X-lens: a message 'Change immersion" appears on the screen. Put a **little bit** oil on your slide, press 'done' and focus again



(IEN) TI	les - Ad	vanced	Setup - Z	IEN 3.7	(ZEN s	ystem			
File	Edit	View	Acquis	sition	Grap	hics	Tools	Window	Help
@	M			Save	All	†m			
رم Loc	ate	Acqui] sition	Proces] sing	Ai	nalysis	Applicati	ons 👻
Transr	nitted	Light	Off			Refle	ted Ligh	nt Off	On
Switch	1 To	Ø	Transn	nission)		6	* Fluores	cence
Favori	tes [Configu	re						
A	I Off	C	API	4	88	1	561		
100	≝ Mici	oscope	Contro	1			3	Show All	

9. Switching to other lenses from an oil-lens or from a dry lens to an oil-lens will result in a message 'Change immersion': please apply immersion or clean objective/specimen'. After applying oil or removing oil press 'Done'.

BE SURE TO REMOVE THE OIL WHEN SWITCHING TO THE 20X-LENS!!



Making a new configuration and a quick first confocal image

Tip: undock the channels window by clicking on the arrow pointing up to the right

- 1. Click 'acquisition' and pull the slider on the microscope totally out (right side of the oculars on the microscope)
- 2. Click 'Smart setup'
- 3. Click on '+Dyes & Contrast Methods'
- 4. Double-click on a dye. The dye will appear in the Smart Setup
- 5. To set up your acquisition channels you have to select all the dyes that are in your sample.
- 6. When done, close the 'Add dye or contrast Method' window
- 7. If you want you can change the colors of the channels by clicking on the arrowhead next to the color
- Choose one of the proposals (fastest, best signal, smartest) and click 'OK'. Best signal takes longer but gives better results (less

crosstalk). Smartest (line-switching) is fast but also very accurate.

- 9. Go to the 'Acquisition Mode'-window, choose an 'Averaging number'. The higher the number, the less noise but also the longer it will take to make the image: quality vs. time, your decision.
- 10. Click on the button 'Confocal': sampling will be set to 1 and the frame size will be adjusted to get the best pixel size calculated from the numerical aperture and Nyquist rate. You can always change the frame size by clicking on 'Presets' and choose another setting.
- 11. Go to the channels window and turn one track on for focussing
- 12. Click 'Set exposure'
- 13. Click 'Live' and adjust the focussing
- 14. Turn on all tracks and click 'Set exposure'
- 15. Click 'Snap': a first confocal image will appear.



Automated Image Export Experiment Feedback Bleaching Direct Processing ZEN Data Storage Upload Automation Acquisition Mode Show All Acquisition Mode Show All Image Size 319.5 µm O Scan Area IDX Image Size 319.5 µm O Scan Area IDX Image Size 319.5 µm Prame Inte Frame Size 2586 px Sampling 1.0 x O Scan Area Image Size Image Size 319.5 µm Prame Time 51.16 s Pixel Time 0.41 µs Scan Speed Image Size Image Size 1.0 000° Correction X 0.000° Line Step Image Size Averaging None Wode Repeat per Line Repeat per Line Repeat per Frame Mode HDR Illumination Wolds HDR Illumination	Experiment R	egions 🔲 Auto Save	Histo
I Bleaching Direct Processing ZEN Data Storage Upload Automation Acquisition Mode Show Al 2 LSM Frame Line Crop Area 0 Scan Area Image Size 319.5 µm × 319.5 µm Pixel Size 0.12 µm Frame Size 2586 px : × 2586 px : Presets Sampling 1.0 x Confocal Frame Time S1.16 5 Pixel Time 0.41 µs Scan Speed T Max Direction Auto Correction X 000 C Correction X 000 C Correction X 000 C Line Step Averaging None 2x 4x 8x 16x Mode Repeat per Line Repeat per Fram. Method Mean Intensity Sum Intensity Bits per Pixel 8 16 Channel HDR Illumination	Automated In	nage Export 🛛 Experiment Feedback	
ZEN Data Storage Upload Automation	Bleaching	Direct Processing	Coloca
Acquisition Mode Show All Acquisition Mode Show All	🗌 ZEN Data Sto	rage Upload 🛛 Automation	#
Acquisition Mode ✓Show All ⊄ LSM Frame Line Crop Area O Scan Area Image Size 319.5 µm 9ixel Size 0.12 µm Frame Size 2586 px ÷ × 2586 px ÷ Presets * Sampling 1.0 × Confoca Frame Time 51.16 5 Pixel Time 0.41 µs Scan Speed			Unmb
ISM Frame Line Crop Area 0 Iox : Ix • Scan Area Iox : Ix Iox : Ix Image Size 319.5 µm × 319.5 µm Pixel Size 0.12 µm Frame Size 2586 px : × 2586 px : Presets * Sampling 1.0 x Confoca Frame Time 51.16 5 Pixel Time 0.41 µs Scan Speed Imax Imax Direction Image Size Correction X 0.000 * Image Size Zoom Correction X 0.000 * Image Size Commensity Immensity Immensity Node Repeat per Line Repeat per Fra Tools Channel HDR Illumination # Channels Image Size Image Size Image Size Image Size Image Size Image Size	Acquisiti	on Mode Show All 🖉	
Frame Line Frame Line O Scan Area JOX TX Image Size 319.5 µm × 319.5 µm Pixel Size 0.12 µm Frame Size 2586 px T × 2586 px Presets * Sampling J.0 X Confocal Frame Time 51.16 s Pixel Time 0.41 µs Scan Speed T Max 1 Direction Correction Auto Coalog T Max 1 Dimension Averaging None 2x 4x 8x 16x Zoom Mode Repeat per Line Repeat per Fra Mode Channel HDR Illumination Averaging HDR Illumination Image Size Image Size Image Size Image Size	ISM		
Frame Line Crop Area Q O Scan Area Image Size 319.5 µm Image Size 319.5 µm Prame Size 2586 px 2586 px > Frame Size 2586 px Sampling 1.0 x Confocal Frame Time 51.16 s Pixel Time 0.41 µs Scan Speed 7 Max Image Size Direction Image Size Correction X 0.000 Correction X 0.000 Correction X 0.000 Correction X 0.000 Mode Repeat per Line Repeat per Line Repeat per Fra Method Mean Intensity Bits per Pixel 8 HDR Illumination Channel	LOIVI	i	i
Crop Area Crop A	Fi	ame Line	
O Scan Area Image Size 319.5 µm × 319.5 µm Pixel Size 0.12 µm Frame Size 2586 px * 2586 px * Presets * Sampling 1.0 x Confoca Frame Time 51.16 5 Pixel Time 0.41 µs Scan Speed	Crop Area	Q (<u>1.0 x</u> ∓ 1.x	
Image Size 319.5 µm × 319.5 µm Pixel Size 0.12 µm Frame Size 2586 px > 2586 px Presets * Sampling 1.0 x Confocal Frame Time 51.16 s Pixel Time 0.41 µs Scan Speed Image Size Image Size Image Size Image Size Direction Image Size Image Size Image Size Image Size Image Size Direction Image Size Image Size Image Size Image Size Image Size Image Size Direction Image Size Image Size Image Size Image Size Image Size Image Size Direction Image Size Image Size Image Size Image Size Image Size Image Size Direction Image Size Image Size Image Size Image Size Image Size Image Size Direction Image Size Image Size <td< td=""><td>O Scan Area</td><td>and the second second</td><td></td></td<>	O Scan Area	and the second	
Image Size 319.5 µm × 319.5 µm Pixel Size 0.12 µm Frame Size 2586 px × 2586 px Presets * Sampling 1.0 x Confoca Frame Time 51.16 5 Pixel Time 0.41 µs Scan Speed			
Frame Size 2586 px > 2586 px Presets * Sampling 1.0 x Confocal Frame Time 51.16 s Pixel Time 0.41 µs Scan Speed 7 Max i Details * Direction 7 Max i Details * Direction Correction X 0.000 ° correction X coord * Averaging None 2x 4x 8x 16x Zoom Mode Repeat per Line Repeat per Fra Tools Channels Channels # Channels # 16 16 Channels Channels Channels	Image Size	319.5 μm × 319.5 μm Pixel Size 0.12 μm	
Sampling 1.0 x Confocal Frame Time 51.16 5 Pixel Time 0.41 µs Scan Speed 7 Max Direction 2 Correction X 0.000° Correction Auto Correction Y 0.000° Line Step 1 Averaging None 2x 4x 8x 16x Correction Y 0.000° Line Step 1 Averaging None 2x 4x 8x 16x Mode Repeat per Line Repeat per Fra Mode Mean Intensity Sum Intensity Bits per Pixel 8 16 HDR Illumination Channels	Frame Size	2586 px + × 2586 px + Presets *	
Frame Time 51.16 s Pixel Time 0.41 µs Scan Speed	Compling	10.	
Frame Time \$1.16 s Pixel Time 0.41 µs Scan Speed 7 Max i Details i Direction Auto 0.000 ** Correction X 0.000 ** i Line Step 1 * Averaging None 2x 4x 6x 16x Mode Repeat per Line Repeat per Fra Tools Channels Bits per Pixel 8 16 Channels Channels A channels 1 1 1 1	Sampling	1.0 X Confocal	
Frame Time 51.16 s Pixel Time 0.41 µs Scan Speed Image: Constraint of the system of the	Late out		
Scan Speed 7 Max Direction 2 Details * Direction Auto Correction X 0.000 C Correction X 0.000 C Correction X 0.000 C Line Step 1 C Averaging None 2x 4x 8x 16x Zoom Mode Repeat per Line Repeat per Fra Mode Mean Intensity Sum Intensity Bits per Pixel 8 16 Channel HDR Illumination	Frame Time	51.16 s Pixel Time 0.41 µs	
Direction Correction X Correction X Correction Y Core	Scan Speed	T Max	
Direction Correction X Correction X Correction X Correction Y Correction X Correcti	1	Details »	
Direction Correction Auto Correction Correction X Correction X Correction V 0.000 Correct	-		
Correction Auto Correction X 0.00° Correction X 0.00° Correction Y 0.	Direction		
Correction X 000° Correction Y	Correction	Auto	
Correction Y 0.00 C Line Step 1 Averaging None 2x 4x 8x 16x Mode Repeat per Line Repeat per Fra Method Mean Intensity Sum Intensity Bits per Pixel 8 16 HDR Illumination Channels	1	Correction X 0.00 ° :	
Line Step 1 Line Step 1 Averaging None 2x 4x 8x 16x Mode Repeat per Line Repeat per Fra Method Mean Intensity Sum Intensity Bits per Pixel 8 HDR Illumination Channels Channels	1	Correction Y 0.00 ° 🗧	
Averaging None 2x 4x 8x 16x Averaging None 2x 4x 8x 16x Mode Repeat per Line Repeat per Fra Zoom Method Mean Intensity Sum Intensity Tools Bits per Pixel 8 16 Channel Image: A channels Image: A channel Image: A channel	Line Sten		
Averaging None 2x 4x 6x 16x Mode Repeat per Line Repeat per Fra Zoom Method Mean Intensity Sum Intensity Tools Bits per Pixel 8 16 Channel HDR Illumination Channels X			Dimensio
Mode Repeat per Line Repeat per Fra Zoom Method Mean Infensity Sum Intensity Tools Bits per Pixel 8 16 Channel HDR Illumination L L	Averaging	None 2x 4x 8x 16x	
Method Mean Intensity Sum Intensity Tools Bits per Pixel 8 16 HDR Illumination Channels	Mode	Repeat per Line Repeat per Fra	Zoom
Bits per Pixel 8 16 Channel Channel Channel	Method	Mean Intensity Sum Intensity	Tools
HDR Illumination	Pite por Divol	16	
HDR Illumination	bits per Pixer		Channel
🔺 Channels 🖬		HDR Illumination	
🕐 Channels 📓	-		
	👍 Channels	Ľ.	





Zooming in

- 1. In the 'Acquisition Mode'-window click on 'Crop Area': a square will appear on screen
- 2. Click on the little cross, hold and move the square over the object of interest
- 3. The square is already zoom 2. You can increase the zoom by clicking on a corner, hold it and move the mouse so a smaller square will be made. Another way is using the slider or arrowheads next to 'Crop Area'
- 4. The frame size will be automatically adjusted, to get the best pixel size.
- 5. Click 'Live' to see if the image is OK or/and click Snap'.

Adjusting tracks manually

Set-exposure can result in too high signals for some specimens. You have to adjust the settings by hand

- 1. Make a channel active in the 'Channels'- window. You can leave all tracks on!
- 2. Click live and check on the 'Range indicator'. A black and white image appears where blue has the grey value 0 and red has the grey value 255 (overexposed)
- 3. When necessary put the 'Digital Gain' back to 1
- 4. When necessary lower the 'Digital Offset' to a negative value to put the background to 0. Normally the current value (0) is OK. It is also possible to get rid of auto-fluorescence or a-specific staining by lowering the 'Digital Offset' more. But don't go too low or you will remove your positive signal!
- 5. Change the 'Master Gain' to a value that your positive signal has just a few red pixels.
- 6. The higher you put the 'Master Gain' the more noise will appear.
- 7. Avoid the horizontal stripes that appear when the 'Master Gain' is set too high! When this happens, raise the 'Digial Gain' to make weak signals visible
- 8. You can also raise the laser-power to get more signal, but don't go too high or you will bleach your slide. This differs per fluorochrome.
- 9. When done, activate another channel and do the adjustments again.
- 10. **Be aware:** when using line switching per track instead of frame-switching, the channels that are shown on the same detector (see the tracks in the 'Imaging Setup'-window) have the same 'Master Gain' setting. Set a common gain for both channels and do a fine adjustment with the laser power.
- 11. When all channels are set, click 'Snap'. When switching to a different lens, check the settings again. They will also change between lenses.

A Channels			Show All	Ľ
Track1 Confe	cal AF555		Ref. 📕	-
Track2 Confe	cal AF488			
Track3 Confe	ocal DAPI			
	AF647			•
✓ ∧ + ₫	j Focus Ref.		*	
🗏 High Intensity L	aser Range			
Track1				
Lasers	405 488	🗹 561 🗌	640	
561 nm		-0	2.5 %	٢
Pinhole	-0		23 µm	•
0.82 Airy Units ≙ 1	.1 μm section		1 AU Ma	x
Alexa Fluor 555	-	_		
Master Gain		0	918 V	•
Digital Offset		0-	0	•
Digital Gain	-0		1.0	•
Display Setting	Default			



Saving the image

There are 2 ways of saving your image in .CZI format: uncompressed or compressed

Uncompressed:

- 1. Click 'File' and choose 'Save as CZI'
- 2. Go to the right network-folder, give your image a name and press 'Save' Compressed:
- 1. Go to 'File' and choose 'Save with options'
- 2. Go to the right network-folder, give your image a name
- 3. Choose a compression method at 'Compression' and press 'Save'

Making images with old or the same settings

- Open a CZI-image that was made earlier; NOT a tif or jpg image!
- Go to the left top of the screen next to 'Smart setup' and press 'Reuse'.
- Activate a track, press 'Live' and if necessary adjust 'Master Gain', 'Digital Offset' and if necessary 'Digital Gain' per Track. Press 'Stop' and go to the next track for adjusting.
- Check on all tracks and press 'Snap'

Making a Z-stack

- 1. Check in 'configuration' the 'Z-stack' on
- 2. Click on the 'Z-stack' bar: the settings appear
- 3. At 'Interval' set a value; this is the space between 2 consecutive slides. If you press 'Optimal' that value will be taken over in 'Interval' (half of the pinhole-setting)
- 4. Press 'Center', the other 'Center', press "First last'
- 5. Activate the track with the biggest volume
- 6. Click 'Live'
- Move the fine coarse adjustment knob of the microscope forwards until there is no more signal. The figure in 'Zstack' shows a blue square going up towards 'L'.
- 8. Click 'Set last'
- 9. Move the fine coarse adjustment knob of the microscope towards you until there is no more signal. The figure in 'Z-stack' shows a blue square going down towards 'F'.
- 10. Click 'Set first'; the number of slices will be automatically set
- 11. Move the fine coarse adjustment back to put the current level to the middle and click 'Stop'
- 12. Select "full Z-Stack per Track" to speed up the acquisition
- 13. Be sure all tracks are on, set 'Averaging' to 2 or 4 and click 'Start experiment'
- 14. A series of images will be taken with a distance of 'interval' between every slice. Save the series. On the left of the image a series of buttons appear (2d, Gallery, 3D etc.): try them out, see what happens. Nothing can go wrong.
- 15. In 'Dimensions' a slider appears at 'Z-position for stepping through the series



Z-Stack	43 Slices
🗹 Tiles	4 Tiles
🗌 Time Series	
Full Z-Stack pe	r Track 🔹



Tile module

See Picture A next to microscope for yellow numbers

Making an overview: Decide if you want to make an overview. If not: go to Tile-scan with high magnification

- 1. Decide if you want to make an overview with the 5X-lens (mostly used to see where tissue or cells are with 1 color) or the 10X-lens (more detailed information and better if you want to see more than 1 color)
- 2. Change to the correct lens (make sure you remove any oil if necessary)
- Click on the arrowhead next to 'Default-EMC' (1) and open either '5X-overview-diag-coupe' (whole slide, rectangle coverglass), '5X-overview-cells' (round coverglass) or '10X-overview-line-4colors'
- 4. Turn off the colours you don't want to use in 'Channels' (2)
- 5. When necessary check on 'Tiles' (3), open the 'Tiles' window and click on 'Show viewer' (4)
- 6. Zoom out until you see the contours of the slide; the blue cross is where the image will be made at this moment.
- 7. You can move the table by double clicking in the contour of the slide

5X-overview:

- 1. When necessary click on the tile and adjust the size with the white edges (5). You can move the tile by clicking on the cross (6), hold it and move it.
- 2. In 'Channels' (2) activate a track you want to record, press 'Live', focus and adjust the settings. Do this for all tracks you want to record.
- 3. In the 'Tiles'-window right-click on the tile (TR1) (7) and click 'Set current Z for selected tile regions'.
- 4. In the Tile-viewer click 'Preview' (8), check off 'Use experiment settings' (9) and choose if you want to 'Delete Existing Preview Images' (10)
- Turn off the channels (11) you don't want to record by clicking on the name and click 'Start' (12). You can stop the scanning whenever you want.

10x-overview:

- 6. In 'Channels' (2) activate the channel you want to record, press 'Live', focus and adjust the settings.
- 7. Make a new tile by doing one of these options:
- 8. Click on the square with the plus (13) and define a tile by numbers at x and y, click 'Add tile region'
- 9. Click on the square with the plus (13), click on the circle and define a tile by either number of tiles or diameter in um, click 'Add Tile Region'
- 10. Click on the square without the plus (14), choose rectangle, circle or freehand and draw a region
- 11. You can define more than 1 region
- 12. In the tile-window at 'tile-regions' (4) a new tile appears. When the tile is checked off the tile will not be visible in the tile-viewer and will not be recorded.
- 13. In the 'Tiles'-window right-click on the tile (TR+nr, 7) and click 'Set current Z for selected tile regions'. When you have more regions/tiles you can click "verify" and check the Z-settings for every region or tile.
- 14. In the Tile-viewer click 'Preview' (8), check off 'Use experiment settings' (9) and choose if you want to 'Delete Existing Preview Images' (10)
- 15. Turn off the channels (11) you do not want to record by clicking on the name and click 'Start'.

7



 Images and Documents
 [2]

 Container 1
 Tiles ...d Setup

 - / Image=-05*

 - / Image--05*

 - / Image--06*

 - / Image--06*

 - / DAzeis_ive.czi

 256 kB / 6.79 MB
 01-u_sed.zzi

 01-u_sed.zzi
 C ≈

 - Z21.42 MB / 223.41 MB
 E

When you check off the tile in the Tiles-window (7) you will only see the preview image. When you double click on an area in that image the stage will go to that area.

If you want to make a tile-scan with a higher magnification:

Tile-scan with high magnification

- 1. Change the lens
- 2. Make a new configuration (smart setup) or load an image and press 'Reuse' (right below 1)
- 3. If necessary change the image settings in the 'Acquisition Mode' window
- 4. Check on 'Tiles' (3) and click 'Show viewer' (4) in the Tiles-window
- 5. If necessary right-click in the viewer, choose 'Import preview image...' and load your previously saved preview image
- 6. Double-click on an area you want to image or go to an area by eye (locate)
- 7. Click 'Live' and adjust the focus and settings (laser, 'Master Gain' etc.) of all channels you want to record
- 8. Make a new tile by doing one of these options:
- 9. Click on the square with the plus (13) and define a tile by numbers at x and y, click 'Add tile region'
 - a. Click on the square with the plus (13), click on the circle and define a tile by either number of tiles or diameter in mm, click 'Add Tile Region'
 - b. Click on the square without the plus (14), choose rectangle, circle or freehand and draw a region; right-click for ending freehand drawing
 - c. In the 'Tiles'-window right-click on the tile (TR+nr, 7) and click 'Set current Z for selected tile regions'. When you have more regions/tiles you can click "verify" and check the Z-settings for every region or tile. When you have very big images which are not completely straight, you can add multiple focus points (see below).
- 10. You can define more than 1 region
- 11. Check on 'Automation' (15)
- In the automation-window check on 'Run macro after experiment execution', click on the dots and choose '01 Automated Stitching Active-Image.py'
- 13. Start the imaging by clicking 'Start Experiment' (16).
- 14. Save the result as CZI

🗧 Au	tomation	Ċ
🔳 Run r	nacro before experiment execution	
Macro		
🗹 Run r	nacro after experiment execution	
Macro	01 Automated Stitching Active Image.pv	



The result images can also be reloaded in the tile-viewer in a later stage, even on another day. Be sure to put the right slide in the same way under the microscope!

Images with more scenes can be split into separate images at 'Processing'-'Utilities'-'Split Scenes', also possible with the freeware Zen-software you can install on your own computer.

Adding multiple focus points

- 1. In the tile module click "Focus Surface and Support Points"
- 2. Make the tile where you want to add support points active by clicking on the name
- 3. Scroll down to "Add Multiple Support Points"
- 4. There are 2 ways of adding support points:
 - a. Generic: choose the number of coloumns and rows and press 'Distribute'
 - b. Onion skin: choose density and margins and press 'Distribute'.
- 5. When you are not happy with the distribution you can press

is the second se

- 6. In the Tile small circles with dots appear that indicate the focus points.
- 7. You can verify the support points or do an autofocus before scanning

Verify

- 8. Click "Verify" and click "Live" to see the live image.
- 9. Click "Move to Current Point" and set the correct focus. When the focus of this point is correct, press "Set Z & Move to Next". The system will go through all the points and remember the Z-position.
- 10. When all points have been verified, click close and press Stop.

Autofocus:

- 11. Open the 'Focus strategy' window
- 12. At the first arrowhead choose 'Use Z values/Focus surface defined in Tiles setup'
- 13. At the second arrowhead choose 'By software autofocus'









- 14. Open the 'Software autofocus' window
- 15. Activate 'Intensity', 'Smart', 'Default', 'Relative Range' and check on 'Automatic Range'

Now you can image as before:

- 16. Check on 'Automation' (15)
- In the automation-window check on 'Run macro after experiment execution', click on the dots and choose '01 Automated_Stitching_Active-Image.py'
- 18. Start the imaging by clicking 'Start Experiment' (16).
- 19. Save the result as CZI

It is possible to combine a tile-scan with z-stacks! Be aware of the time this will take.

Shutting down the system

- 1. On the microscope touchscreen press 'Load position'
- 3. If necessary clean the oil lenses
- 4. Save the images you want to save
- 5. Close the Zen software
- 6. Have a look in the agenda and see if there will be someone after you
- 7. Someone after you:
- 8. Leave the system as it is
- 9. Last user of the day:
- 10. Shut down the computer
- 11. Turn off the connector box on the microscope table
- 12. Turn off the 'system'-switch behind the table, box above the box with the key. DO NOT TOUCH OTHER SWITCHES OR THE KEY!



- Au	tomation	1
💷 Run n	nacro before experiment execution	
Macro		